QUANTOFIX® Kalium

de

Inhalt:

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en

Pack content:

- 1 aluminum container with 100 test strips
- 2 bottles K⁺-1
- 2 test tubes

Measuring range:

200 – 1500 mg/L

Color gradation:

0 - 200 - 400 - 700 - 1000 - 1500 mg/L

General indications:

Remove only as many test strips as are required. Close the container immediately after removing a strip. Do not touch the test field.

Additional indications:

The orange-red coloration of the unused test fields is normal and does not indicate any decomposition of the reagents.

Instructions for use:

- Place the test tube into the cavity of the thermoformed mold.
 - Add 10 drops K⁺-1 (0.7 % nitric acid).
 - Dip the test strip into the neutral sample for 1 second.
 - Shake off excess liquid.
 - Place the test strip into the test tube.
 - Wait 60 seconds.
 - Shake off excess liquid.
 - Compare test field with the color scale.
- If potassium ions are present, the test field turns yellow to orange-red.

Interferences:

Sodium ions in concentrations > 3000 mg/L lead to a fading of the color reaction with corresponding low test results. The following ions in the indicated concentrations do not interfere:

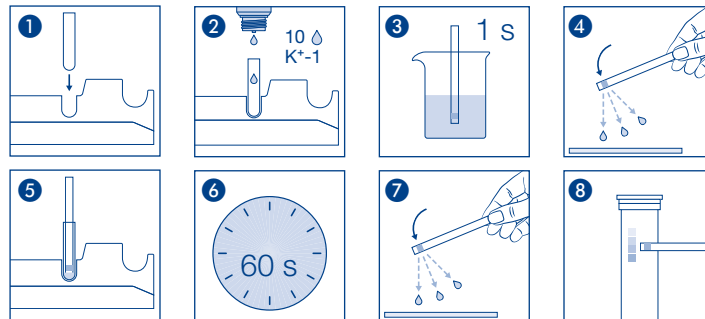
- < 1000 mg/L Al³⁺, Ba²⁺, Bi³⁺, Ca²⁺, Fe^{2+/3+}, Mg²⁺, Mn²⁺, Sr²⁺, Zn²⁺
- < 200 mg/L NH₄⁺, Hg²⁺, Ti⁺
- < 25 mg/L S²⁻

Disposal:

Information regarding disposal can be found in the safety data sheet. You can download the SDS from www.mn-net.com/SDS.

Storage:

Avoid exposing the strips to sunlight and moisture. Keep container cool and dry (storage temperature 4 – 30 °C). If correctly stored, the test kit may be used until the use-by-date printed on the packaging.

MACHEREY-NAGEL GmbH & Co. KG · Valenciener Str. 11 · 52355 Düren · Germany
Tel.: +49 24 21 969-0 · info@mn-net.com · www.mn-net.com

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MACHEREY-NAGEL GmbH & Co. KG · Valencienner Str. 11 · 52355 Düren · Germania